Microscope · Mikroskop · Microscope

DST-1028

**EN** Instruction manual

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**EN** Visit our website via the following QR Code or web link to find further information on this product or available translations of these instructions.

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**DOWNLOAD "CamLabLite" SOFTWARE FOR WINDOWS**

[archive.bresser.de/download/CamLabLite/](http://archive.bresser.de/download/CamLabLite/)

<b>English</b> .....	<b>4</b>
<b>Deutsch</b> .....	<b>10</b>
<b>Français</b> .....	<b>16</b>

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For any warranty claims or service inquiries, please refer to the information on "Warranty" and "Service" in this documentation. We ask for your understanding that unsolicited returns cannot be processed.

Errors and technical changes excepted.

© 2022 Bresser GmbH

All rights reserved.

The reproduction of this documentation - even in extracts - in any form (e.g. photocopy, print, etc.) as well as the use and distribution by means of electronic systems (e.g. image file, website, etc.) without the prior written permission of the manufacturer is prohibited.

The designations and brand names of the respective companies used in this documentation are generally protected by trade, trademark and/or patent law in Germany, the European Union and/or other countries.

## 2 Validity note

This documentation is valid for the products with the following article numbers:

5808600

**Manual version:** 0522

**Manual designation:**

Manual\_5808600\_DST-1028-USB-Microscope\_en-de-fr\_BRESSER\_v052022a

Always provide information when requesting service.

## 3 About this Instruction Manual

### NOTICE

**These operating instructions are to be considered a component of the device.**

Read the safety instructions and the operating manual carefully before using this device.

Keep this instruction manual in a safe place for future reference. When the device is sold or given to someone else, the instruction manual must be provided to the new owner/user of the product.

---

## 4 Intended use

- This device is not intended for use by persons (including children) with impaired physical, sensory or mental abilities, or for lack of experience and/or knowledge, unless they are supervised by or have received instructions from a person responsible for their safety on how to use this device.
- This device is for personal use only.
- It was developed for an magnified representation of nature observations.
- The device is intended only for indoor use.

---

## 5 General safety instructions

### DANGER

#### **Danger of suffocation!**

Improper use of this product may result in suffocation, especially for children. It is therefore imperative that you observe the following safety information.

- Keep packaging materials (plastic bags, rubber bands, etc.) away from children! There is a danger of choking!
- This product contains small parts that can be swallowed by children! Choking hazard!
- If small parts are swallowed, consult a doctor immediately!

### CAUTION

#### **Risk of body injury!**

This device contains components and/or accessories that can cause minor or serious body injury if used improperly. It is therefore essential that you observe the following safety information in order to prevent body injuries.

- Sharp-edged and pointed tools are often used for working with this device. Therefore, keep this device and all accessories and tools in a place inaccessible to children. Danger of body injury!
- Chemicals and liquids do not belong in children's hands! Don't drink! Clean hands thoroughly under running water after use. In case of accidental contact with eyes or mouth, rinse with water. In case of complaints, consult a doctor immediately and show the substances.

### **NOTICE**

#### **Danger of material damage!**

Improper handling may result in damage to the unit and/or accessories. Therefore, use the device only in accordance with the following safety information.

- Do not disassemble the device! In the event of a defect, please contact your dealer. They will contact the Service Center and can arrange the return of this device for repair if necessary.
- Do not expose the device to high temperatures and protect it from water and high humidity.
- Do not subject the device to excessive vibrations.
- Only use accessories and spare parts for this device that comply with the technical specifications.

## 6 About this microscope

This microscope is only a digital microscope and therefore does not offer the possibility of visual observation.

It is equipped with an objective lens and a camera sensor for image recording. The digital images can be transferred directly to a Windows PC via the supplied USB cable, being viewed there and evaluated, optimised and finally saved using the supplied software\*.

\*Windows version available for download (see download info on p. 2)

## 7 Choice of location

Choose a suitable location before installation and commissioning.

Place the device on a stable, level and vibration-free surface.

## 8 Parts overview and scope of delivery

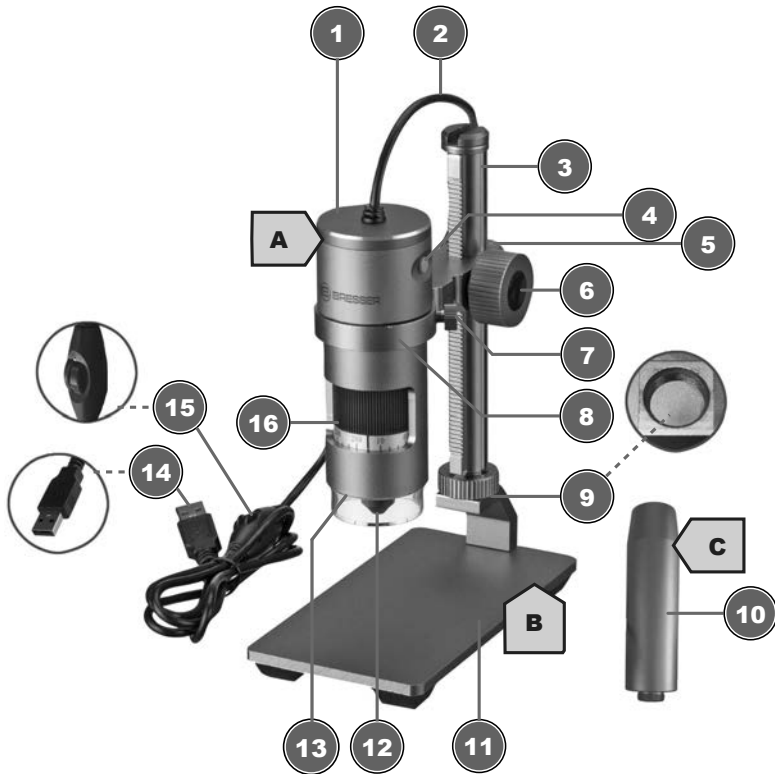
Illustration 1: All parts of the digital microscope DST-1028.

1	Microscope head	2	USB connecting cable
3	Running rail with integrated rack	4	Trigger <sup>1</sup>
5	Locking screw (holder)	6	Wheel for height adjustment / focusing
7	Fixing screw (holder)	8	Digital microscope holder
9	Holding thread	10	Extension piece
11	Microscope base	12	Objective
13	LED ring lighting	14	USB-A connector
15	Dimmer wheel	16	Zoom ring with scale (10x to 280x)

<sup>1</sup>Function not available for this model

### Scope of delivery:

Digital microscope (A) with USB cable, microscope stand with running rail and holder (B), extension piece (C)

---

## 9 Assembling the microscope

This microscope is designed for operation with a Windows PC.

For the assembly proceed as follows:

1. Screw the running rail with the preassembled microscope holder into the holding thread. Make sure that the holder is facing forward as shown.
2. **NOTICE! For higher observation objects, the supplied extension piece can first be screwed into the mounting thread. The rail is then screwed into the extension piece. This will increase the working height significantly.**
3. Loosen the fixing screw on the holder and unscrew it slightly so that it does not block the retaining ring.
4. Insert the microscope head into the holder and hand-tighten the fixing screw.

## 10 Power supply

Plug the USB-A connector of the USB cable into a free USB port of the PC. The power is supplied through the computer after it has been switched on. The LED ring lighting will light up.

## 11 Connecting to a PC

Use of this digital microscope requires a connection via USB to a PC running the Windows 10 operating system. Follow the steps below for proper connection:

1. Download and install the CamLabLite\* software on the PC.
2. Mount the digital camera as described.
3. Insert the USB-A plug of the USB connection cable into a free USB port of the PC. If the connection is successful, the LED lights up. The digital microscope is ready for use.
4. The CamLabLite\* software can be used to view and save images from the digital microscope. Furthermore, various image optimizations can also be made. Please also read the separate software instructions.

*\*Information about downloading the software and the software instructions can be found on page 2 of this manual.*

## 12 Use of the software

For this digital microscope, we offer extensive software for editing and saving your digital images. This can be downloaded from the product page via the following link ("Software" section).

Furthermore, you can also find detailed software instructions in different languages under this link (section "Instructions").

<https://www.bresser.de/P5808600>

### **System requirements:**

PC/notebook with Intel i3, i5, i7 or i9 processor from 2nd generation or AMD FX or Ryzen processor, operating system Windows 8.1/10/11, free USB 2.0 port, internet connection for software download

## 13 Adjusting the lighting

The digital microscope is equipped with a LED ring lighting for optimal object illumination and can be dimmed continuously.

1. Connect the digital microscope to the PC.
2. Once the LED lighting is on, the brightness can be adjusted continuously by turning the dimmer with on the USB cable.

---

## 14 Adjusting the height of the head

Depending on the size of the observation object or when changing the magnification setting, the position of the microscope head must be adjusted in height. To do this, proceed as follows.

1. Hold the microscope head and loosen the locking screw for the holder.
2. Turn the height adjustment wheel to move the microscope head up or down in the holder.
3. Hand-tighten the locking screw when the desired height position is reached.

**NOTICE! Especially when changing the zoom range, it may be necessary to additionally change the position of the digital microscope in the holder. (see also "Focusing the object")**

## 15 Adjusting the zoom

The microscope optics have a zoom range from 10x to 280x.

The zoom range can be changed by turning the zoom ring. The corresponding digit on the scale must be visible above the marking arrow.

**NOTICE! When changing the zoom range, the focus may also need to be adjusted. (see also "Focusing the object")**

## 16 Focusing the object

For optimal image sharpness, the observation object must be brought into the focus range of the digital camera. To do this, proceed as follows:

1. Depending on the object height, assemble the extension piece if necessary.
2. Set the required head height for the digital microscope.
3. Make the desired magnification setting.
4. If necessary, correct the height and the position of the digital microscope in the holder slightly until the image on the monitor appears sharp.
5. **WARNING! Always hold the digital microscope firmly when loosening the locking screw or the fixing screw on the holder to prevent the digital microscope from sliding down. The digital microscope could hit the microscope base and be irreparably damaged!**

## 17 Technical data

Zoom range:	10x to 280x
Image field:	27 x 48mm (10x) or 0.9 x 1.7mm (280x)
Working distance:	10mm to 110mm
Max. object height:	approx. 97mm (10x) to 160mm (280x)
Illumination:	8 LEDs
Digital camera:	built-in (Sony IMX335 CMOS sensor with 5.1 megapixels)
PC connection:	USB-A*
Power supply:	5V via USB connection*
<b>Materials</b>	
Objective lens:	glass, multi-coated
Tripod	metal
*Data transfer and power supply are carried out via the USB port on the PC at the same time	

---

## 18 EC declaration of conformity

A "Declaration of conformity" in accordance with the applicable directives and corresponding standards has been prepared by Bresser GmbH. The full text of the EC declaration of conformity is available at the following Internet address: [www.bresser.de/download/5808600/CE/5808600\\_CE.pdf](http://www.bresser.de/download/5808600/CE/5808600_CE.pdf)

## 19 UKCA Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the UKCA declaration of conformity is available at the following internet address: [www.bresser.de/download/5808600/UKCA/5808600\\_UKCA.pdf](http://www.bresser.de/download/5808600/UKCA/5808600_UKCA.pdf)

**Bresser UK Ltd.** • Suite 3G, Eden House, Enterprise Way, Edenbridge, Kent TN8 6HF, Great Britain

## 20 Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!  
According to the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment and its transposition into national law, used electrical equipment must be collected separately and recycled in an environmentally sound manner.

## 21 Warranty

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at [www.bresser.de/warranty\\_terms](http://www.bresser.de/warranty_terms).

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**GB IE**

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

**e-mail:** [service@bresseruk.com](mailto:service@bresseruk.com)  
**Telephone\*:** +44 1342 837 098

BRESSER UK Ltd  
Customer Support  
Suite 3G, Eden House  
Enterprise Way  
Edenbridge, Kent TN8 6HF  
United Kingdom

\*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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